



**Fisherbrand**<sup>®</sup>  
**QUALITY. RELIABILITY. VALUE.**

Versa-Clean<sup>™</sup>

Liquid Concentrate

Readily Biodegradable (OECD 301D).  
 Outstanding dispersing and rinsing properties.  
 Passes Inhibitory Residue Test (Standard Methods for the Examination of Water and Wastewater 20th Ed., 9020B-4A-2).  
 USDA approved for use in Federally inspected meat and poultry plants.  
 Trace Analysis available.

## Environmentally Safe

Does not contain: Alcohol, Ammonia, APE's, Butyl, Chlorine, Caustic, Fragrance, NPE's, NTA, Phosphates, Silicates, Solvents, Sulfates or VOC's, Ozone Friendly.

## User Friendly

Non-Abrasive • Mild pH • Non-Flammable • Non-Fuming  
 Non-Carcinogenic • Pleasant Odor • Streak-Free  
 Safe on Skin • Non-Toxic • Economical Concentrate

## Versatile

Safe for all washable surfaces, harmless to most types of fabrics and metals.  
 Recommended for: aluminum, appliances, chrome, computer screens, countertops, floors, glassware, labware, plastic, porcelain, steel, walls, and much more.

Size	Cat.No.
1L (33 7/8 oz.)	04-342
4L (1 1/16 gal.)	04-343



In Accordance With  
**OECD 301D**



NOT TESTED ON ANIMALS



FULLY RECYCLABLE  
 CARDBOARD AND PLASTIC

SEE REVERSE SIDE FOR USAGE DETAILS  
**BN0317092**

For over 35 years, Fisherbrand Versa-Clean has been used safely and effectively for a wide variety of applications in a host of industries, from laboratories to industrial plants. Eliminates the hazardous concerns associated with petroleum distillates, chlorinated solvents, as well as alkaline and acidic products. Highly recommended for critical cleaning. This emulsifier system incorporates a fine blend of high quality synthetic penetrants and wetting agents.

Aluminum	N/D → 15:1
Appliances	10:1 → 30:1
Blood	N/D → 10:1
Carbon Deposits	N/D → 10:1
Carpets	50:1
Carpets, Spot Cleaning	5:1 → 10:1
Chrome	30:1
Cleaning Equipment	10:1
Computer & Keyboard	30:1
Computer Screens	60:1
Concrete	N/D → 5:1
Diesel Fuel	10:1
Dip Tanks	5:1 → 15:1
Electrostatic Filters	N/D
Floors	40:1
Garbage Containers	N/D → 10:1
Glassware	60:1
Grease Filters	N/D → 10:1

Laundry	1/4 Cup/Load
Laundry Pre-Soak	30:1
Laundry Spots	N/D
Lubricants	N/D → 10:1
Marble	40:1
Oil Deposits	N/D → 10:1
Porcelain	10:1
Pressure Sprayers	30:1 → 50:1
Sinks	N/D → 10:1
Smoke	10:1
Soak Baths	N/D → 60:1
Stainless Steel	30:1 → 50:1
Steam Cleaners	50:1
Steel	N/D → 5:1
Ultrasonic Cleaning	10:1 → 40:1
Vents	N/D → 20:1
Walls	30:1 → 60:1
Wood	30:1 → 50:1

**N/D = NON DILUTED**

The dilution rates serve as guidelines to common applications, in some cases the dilution of water to Versa-Clean may be increased or decreased, according to the severity of the specific application.

## Use With

Automatic Scrubbers • High-Pressure Sprayers  
 Laundry Machines • Metering Equipment • Soak Tanks  
 Spray-on/Wipe-off • Ultra Sonic Cleaners  
 Automatic Glassware and Labware Cleaners

## In a Variety of Industries

Aerospace • Automotive • Chemical Processing • Education  
 Food Processing • Healthcare • Metalworks • Military  
 Research Laboratories



For customer service: call 1-800-766-7000  
 To fax an order: use 1-800-926-1166  
 To order online: [www.fishersci.com](http://www.fishersci.com)

BN0317092